


UVC 500 Ultraviolet Crosslinker

transilluminator

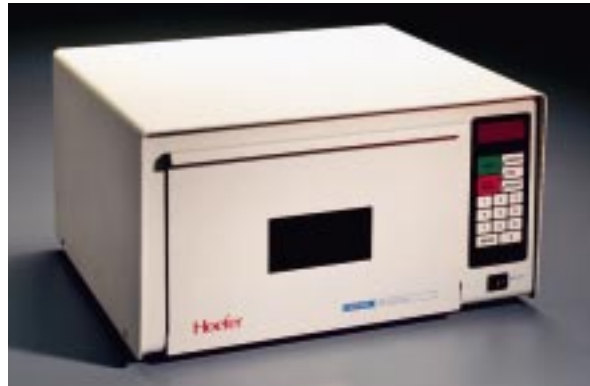

Page finder

UV Crosslinker function and description	1
Unpacking	2
Specifications	2
Important safety information	3
Operating instructions	3
Energy setting	4
Time setting	5
Care and maintenance	6
Bibliography	8
Customer service information	9


UV Crosslinker function and description

The UVC 500 Ultraviolet Crosslinker provides a uniform and intense source of ultraviolet (UV) light for such uses as:

- DNA crosslinking
- DNA nicking
- Rapid Site mapping
- Testing RecA function
- Ultraviolet sterilization
- Elimination of PCR contamination

The internal photo-feedback system continually measures the energy within the chamber and automatically deactivates the UV light source once the programmed energy level is reached. This system also automatically adjusts to variations in UV intensity that occur as lamps age.

Crosslinker front view

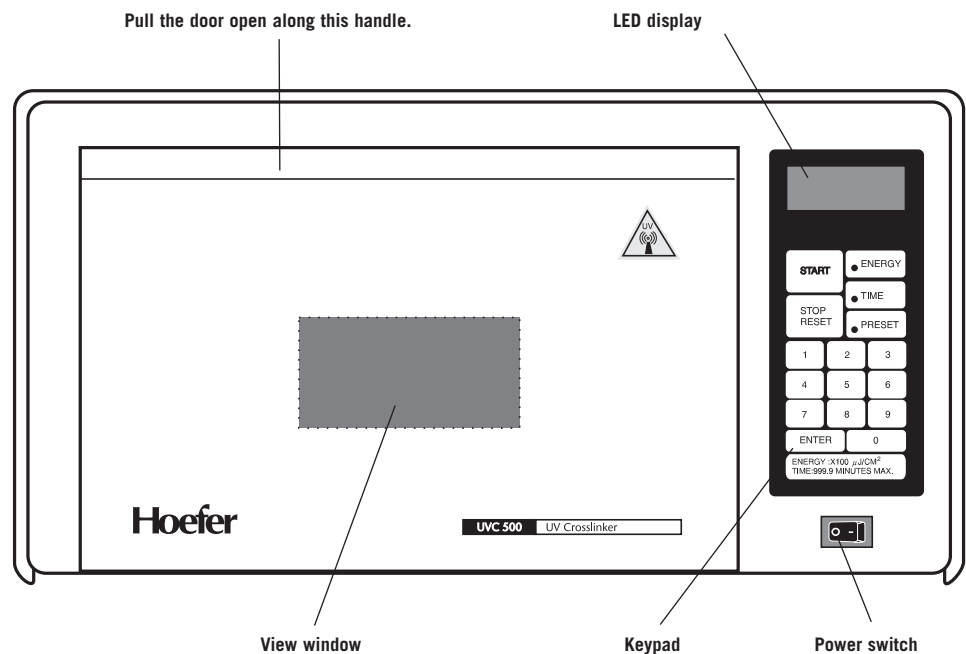

Unpacking

Unwrap all packages carefully and compare contents with the packing list, making sure all items arrived. If any part is missing, contact your local Amersham Biosciences sales office. Inspect all components for damage that may have occurred while the unit was in transit. If any part appears damaged, contact the carrier immediately. Be sure to keep all packing material for damage claims or to use should it become necessary to return the unit.

Specifications

Ultraviolet wavelength and sources

- 254 nm (shortwave)
- Five 8-W 254 nm UV dual bi-pin discharge type tubes

Features

- Microprocessor controlled UV sensor feedback system
- Large LED readout
- Tactile membrane switch keypad
- Laydown-type door
- Multiple set functions: pre-set UV energy or time exposure
user-specified UV energy or time exposure
- Max. energy setting: 999,900 $\mu\text{J}/\text{cm}^2$
- Max. time setting: 999.9 min

Dimensions (h × d × w)

External	22.2 × 34.9 × 40.0 cm (8.75 × 13.75 × 15.75 in.)
Chamber interior	12.7 × 30.5 × 25.4 cm (5 × 12 × 10 in.)

Input fuses

- 115 V - F 1A/250V 5 × 20 mm (2 fuses)
- 230 V - F 1A/250V 5 × 20 mm (2 fuses)


Weight

- 7.5 kg (15.5 lbs)

Safety

- Safety interlocked door
- UV-blocking view window
- Safety-fused and removable power cord

Important safety information

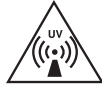

- Always wear protective gloves and goggles when working with the crosslinker.
- The UVC 500 is a powerful source of ultraviolet radiation. The door is safety interlocked to prevent accidental user exposure to UV light. Do not attempt to override the safety interlocks.

If the safety interlock is disengaged, the UV lamps do not light and the unit will beep 3 times when the START key is pressed.

If a malfunction occurs and the lamps in the crosslinker remain on when the door is open, immediately close the door and turn off the mains power. For repair information, call the Amersham Biosciences Technical Support.

- The view window blocks UV radiation and allows viewing while the unit is in operation. If the window is broken or removed, stop using the crosslinker immediately. For repair information, call Amersham Biosciences Technical Support.
- Always disconnect the mains power cord before cleaning the unit or replacing fuses or lamps.

Operating instructions

Set up and membrane placement

1. Place the unit on a stable and level surface.
2. Plug the power cord into the receptacle at the back of the unit and into a properly grounded electrical outlet of the proper power rating.
3. Flip the mains power switch located under the keypad to on (–).
Note: The instrument settings are always the values last programmed, and are indicated by the LED display and a lit red indicator lamp.
4. Open the door by pulling the door handle above the window toward you.
Note: Do not obstruct or damage the photofeedback sensor, which is mounted on a small metal tab protruding from the lower right side.
5. Moisten (do not wet completely) a sheet of blotting paper with transfer buffer and place it in the center of the compartment. Place the membrane on the paper with the nucleic acid side facing upward.
6. Set the energy exposure (energy or time) as described in the next section.

Note: To protect the crosslinker from corrosive buffers, lay a piece of plastic wrap on the compartment bottom before placing the blotting paper and sample.

Energy setting

The amount of energy recommended for UV crosslinking of DNA and RNA by covalently binding nucleic acids to Hybond™ membranes after Northern, Southern, slot or dot blotting is 70,000 $\mu\text{J}/\text{cm}^2$. This value is preset at the factory and displays as 700 on the LED display, which only shows 4 digits. As the unit operates, a countdown of the remaining energy to be emitted is displayed on the LED.

Note: Multiply all displayed energy settings by 100 for the actual energy value.

Using the preset energy setting


1. First press the PRESET key and then the ENERGY key. The red indicator lamp next to each of these keys should light and the LED should display 700 (or the value last programmed).
2. Press the green START key. The UV lamps take a few seconds to energize before the countdown value displays. At the end of the exposure, the unit automatically stops and beeps 5 times.

Reprogramming the preset energy setting

1. Press and hold the PRESET key until you hear the second beep. The lamp next to the PRESET key flashes when this key switches to program mode.
2. Press the ENERGY key. Use the numeric keypad to enter the new value. The LED display should show the new value.
3. Press the ENTER key.

UVC 500 Crosslinker user interface

The LED display shows values for **energy** (multiply by 100 for actual value in $\mu\text{J}/\text{cm}^2$) and **time** (in minutes and tenths of a minute.)

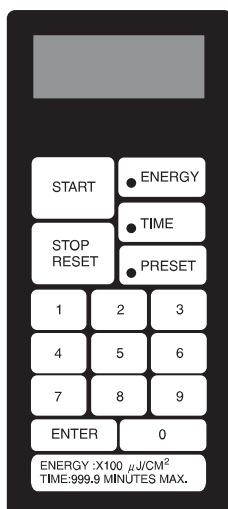

Programming the ENERGY key

1. Press the ENERGY key.
2. Enter a value using the numeric keypad. The LED will display the new value.
3. Press ENTER. After programming, pressing the ENERGY key retrieves this setting. Change the setting by entering a new value and pressing the ENTER key.

Time setting

The display shows time in minutes and tenths of a minute.

Using the preset time setting

1. First press the PRESET key and then the TIME key. The red indicator lamp next to each of these keys should light and the LED should display 5.0 (or the value last programmed).
2. Press the green START key. The UV lamps take a few seconds to energize before the countdown value displays. At the end of the exposure, the unit automatically stops and beeps 5 times.

Reprogramming the preset time setting

1. Press and hold the PRESET key until you hear the second beep. The lamp next to the PRESET key flashes when this key switches to program mode.
2. Press the TIME key. Use the numeric keypad to enter the new value. The LED display shows the new value.
3. Press the ENTER key.

Programming the TIME key

1. Press the TIME key.
2. Enter a value using the numeric keypad. The LED will display the new value.
3. Press the ENTER key. After programming, pressing the TIME key retrieves this setting. Change the setting by entering a new value and pressing ENTER.

Stopping and restarting the crosslinker

- Press the red STOP/RESET key to stop the exposure. The LED will display the remaining exposure level or time.
- Press the green START key to restart the exposure. The program will continue from where it was stopped.
- To enter a different value once the program has started, press STOP/RESET, enter the new value, then press ENTER. Press the START key to begin the new exposure value.
- To revert to the last used setting after the exposure is interrupted, press the STOP/RESET key a second time. Press START to start exposure from the previous setting.
- If the door is accidentally opened during exposure, the UV lamps instantly turn off and the remaining exposure level or time is displayed on the LED display. To restart, close the door and press the START key.

Caution: Always turn the mains power switch off and unplug the instrument before cleaning.

Important: The photofeedback system is calibrated to 254 nm UV light. Do not install lamps rated for a different wavelength.

Care and maintenance

Cleaning

Preventative maintenance includes regular cleaning of the compartment surface, the sensor, and UV lamps.

- Use mild soap and water on a soft cloth or sponge to clean the exterior.
- Gently wipe the photofeedback sensor and UV lamps regularly with a soft cotton cloth dampened with alcohol.
- Always wipe any water and spills from the inside of the unit and clean away all chemicals after each use.
- Never use abrasive cleaners, acetone or chloroform on any part.

Replacing UV lamps

Periodically check that all five lamps function by looking through the view window while the unit is on. If one or more lamps are burned out, it is recommended that all five be replaced at the same time. To replace each lamp:


1. Turn off the mains power switch and disconnect the power cord.
2. Locate the lamps along the top of the compartment. Remove each lamp by carefully rotating it 1/4–1/3 turn and then carefully pulling it loose.
3. To insert the new lamp, hold it so that the metal plugs at the ends align vertically. Place the lamp in the sockets at each end of the unit and lightly push it up into place. Rotate the tube / turn so that it locks into place.

Replacing fuses

Fuses protect equipment by disconnecting loads too large for the instrument's circuit design. For continued protection, only use fuses with the specified voltage and current rating.

Each of two round modules on the left side of the control panel (above the mains power cord receptacle) holds one F 1A/250V 5 × 20 mm (Fast blo) fuse.

UVC 500 Crosslinker back view 115 V model

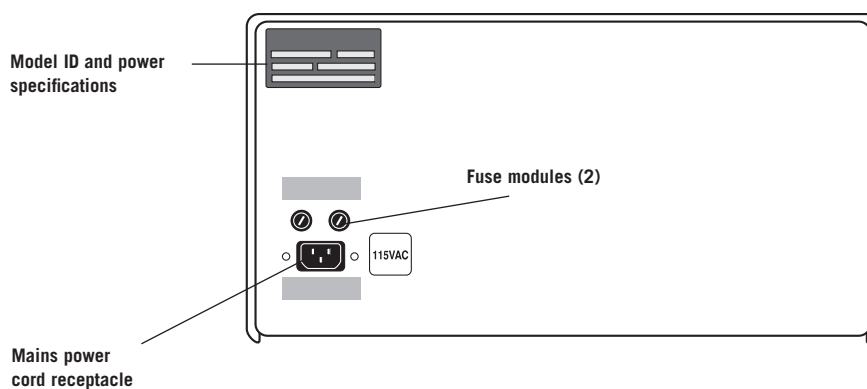

1. **Caution!** Set the mains power switch to off and detach the power cord before replacing fuses.
2. Insert a small flat-blade screwdriver into the slot on the fuse module and turn it 1/4 turn counterclockwise. The spring-loaded module cap will loosen and you can then pull the cap/fuse holder out.
3. Pull the fuse out of its holder and inspect. If the fuse element is burned or broken, replace the fuse with an identical type. If the fuse appears to be intact, check it with an ohmmeter.
4. Insert a good fuse into the holder and then insert this assembly back into the unit. Seat the module by inserting the screwdriver into the slot, pressing gently, and turning the cap 1/4 turn clockwise.

Bibliography

DNA Crosslinking

Church, M. C. and G. Gilbert. (1984) Genomic sequencing. *Proc Natl. Acad. Sci., USA.* **81**, 1991–1995.

Khandjian, E. W. (1987) Optimized hybridization of DNA blotted and fixed to nitrocellulose and nylon membranes. *BioTechnology.* **5**, 165–167.

Khandjian, E. W. (1986) UV crosslinking of RNA to nylon membrane enhances hybridization signals. *Molec. Biol. Report.* **11**, 107–115.

Reed, K. C. and D. A. Mann. (1985) Rapid transfer of DNA from agarose gels to nylon membranes. *Nuc. Acid Res.* **13**, 7207–7221.

DNA Nicking

Vollrath, D. and R. W. Davis. (1987) Resolution of DNA molecules greater than 5 megabases by contour-clamped homogeneous electric fields. *Nuc. Acid Res.* **15**, 7865–7876.

Rapid Site Mapping

Whittaker, P. A. and E. M. Southern. (1986) Ultraviolet irradiation of DNA: A way of generating partial digests for rapid restriction site mapping. *Gene.* **41**, 129–134.

Testing RecA function

Maniatis, T. (1989) Maintenance and Testing of Lysogens of Bacteriophage Lambda in *Molecular Cloning: A Laboratory Manual*, eds. Sambrook, Fritsch and Maniatis. 2.96.

Ultraviolet sterilization

Reich, R. R., and Anderson, H. D. (1986) Sterilization of Membrane Filters with Ultraviolet Radiation. *Pharmaceutical Manufacturing.* Jan.

Elimination of PCR contamination

Wilson, K. H. (1992) UV Absorption Complicates PCR Decontamination, *BioTechniques.* **13**, 208–210.

Sarkar, Gobinda, and Sommer, Steve S. (1991) Parameters Affecting Susceptibility of PCR Contamination to UV Inactivation, *BioTechniques*, **10**, 590–594.

Sarkar, Gobinda, and Sommer, Steve S. (1990) Shedding Light on PCR Contamination. *Nature.* **343**, 27.

Yamada, H. and Hieda, K. (1992) Wavelength Dependence (150-290nm) of the Formation of the Cyclobutane Dimer and the (6-4) Photoproduct of Thymine. *Photochemistry and Photobiology.* **55**, 541–548.

Customer service information

Technical service and repair

Amersham Biosciences offers complete technical support for all our products. If you have any questions about how to use this product, or would like to arrange to repair it, please call or fax your local Amersham Biosciences representative.

Important: Request a copy of the Amersham Biosciences “Contamination Clearance Certificate” before returning the item. No items can be accepted for servicing or return unless this form is properly completed.

Ordering information

	Quantity	Code No.
UVC Ultraviolet Crosslinker		
For membranes to 25 × 30 cm.		
115 VAC	1	80-6222-31
230 VAC	1	80-6222-50
Replacement UV lamp, 8 W, 254 nm	1	80-6223-64
Replacement fuses	5	80-6109-45



Notes:



is a trademark of Amersham Biosciences
Limited or its subsidiaries.

Amersham Amersham Biosciencesand is a trademark of Amersham plc.

All other trademarks and registered trademarks
are the property of their respective companies or
organizations.

© 2000 Amersham Biosciences Inc.
All rights reserved.

All goods and services are sold subject to the terms
and conditions of sale of the company within the
Amersham Biosciences group that supplies
them. A copy of these terms and conditions is
available on request.

Printed in the USA.

Amersham Biosciences UK Limited
Amersham Place Little Chalfont
Buckinghamshire England HP7 9NA

Amersham Biosciences AB
SE-751 84 Uppsala Sweden

Amersham Biosciences Inc.
800 Centennial Avenue PO Box 1327
Piscataway NJ 08855 USA

Amersham Biosciences Europe GmbH
Munzinger Strasse 9 D-79111 Freiburg Germany

www.amershambiosciences.com

